

CEofix™ buffers voor robuste ionen analyse met Capillaire Electrophorese

Analisis R&D s.a.

François de l'Escaille

fde@analisis.be

Jean-Bernard Falmagne

Philippe Louis, Michel Brohet

Studiedag Ionenanalyse KVCV 24 november 2004



analisis

Outline

- Analisis R&D and products
- CEofix™ buffers for CE
- CEofix™ Anions / Cations
- CEofix™ Applications

Analisis R&D laboratory.

For more than 20 years, the R & D (established in Namur) specialized in electrophoresis has developed in vitro diagnosis kits which have been distributed throughout the world.

As a result of our long experience in the development of electrophoresis kits using agarose gels, Analisis has now patented and marketed kits using new automated methods by means of capillary electrophoresis



analisis

Analisis R&D




Bio-clinical Applications

- **Electrophoresis** : ISOPAL, ISOAMYL, PROTUR PLUS, MICROPROTUR, PROTUR B.J., DIATRAC

- **Capillary Electrophoresis**


(P/ACE5000 and P/ACE MDO)


- **HbA1c/HbA2**
hemoglobins and variants
- **SPE/HRE**
serum proteins
- **CDT**
transferrin isoforms (alcohol abuse marker)

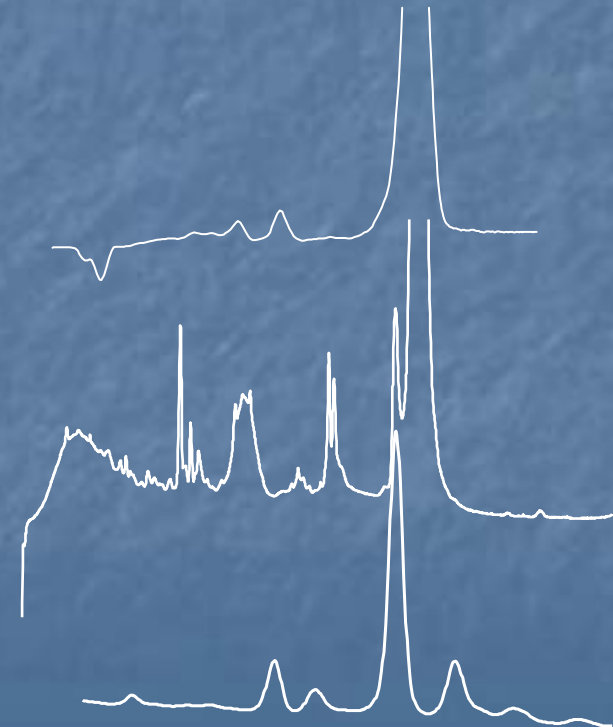

Analytical kits

- Capillary electrophoresis
 - CEofix™ pH
 - CEofix™ MS
 - CEofix™ MEKC
 - CEofix™ Anions
 - CEofix™ Cations
- Editor of PHoEBuS software
(program help for electrophoresis buffer studies)

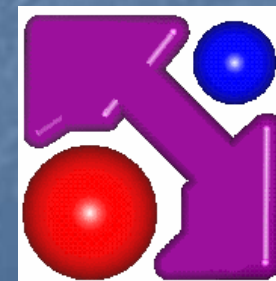


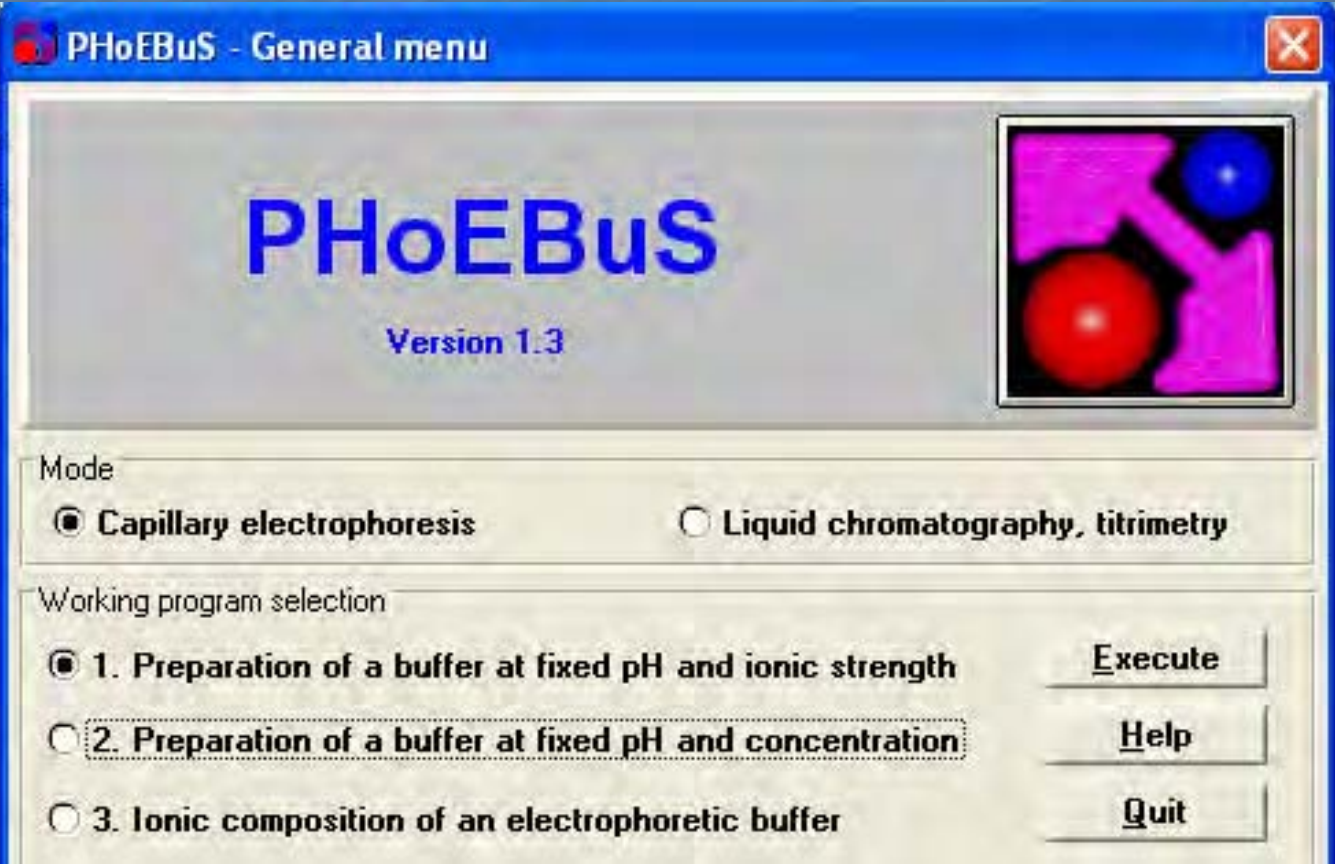

PHoEBuS - General menu

PHoEBuS

Version 1.3

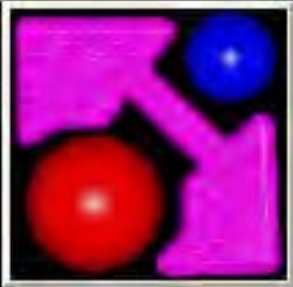

Mode

Capillary electrophoresis

Liquid chromatography, titrimetry

Working program selection

1. Preparation of a buffer at fixed pH and ionic strength

2. Preparation of a buffer at fixed pH and concentration


3. Ionic composition of an electrophoretic buffer

Execute

Help

Quit


Dynamic Coating of capillary

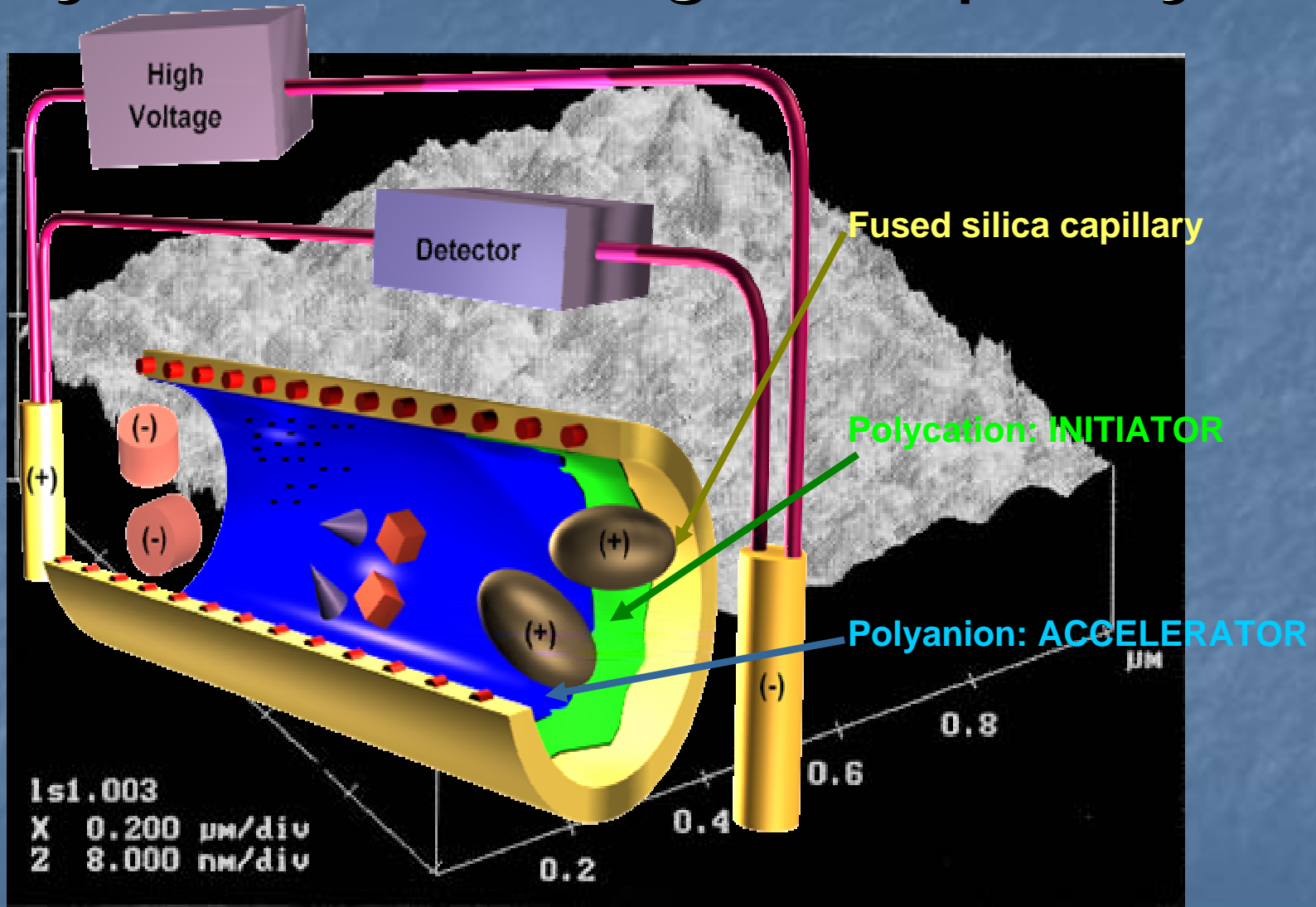

Features of CEofix™

- Control of EOF
- Reproducibility
 - No-memory

Beckman Coulter P/ACE MDQ with 32K software

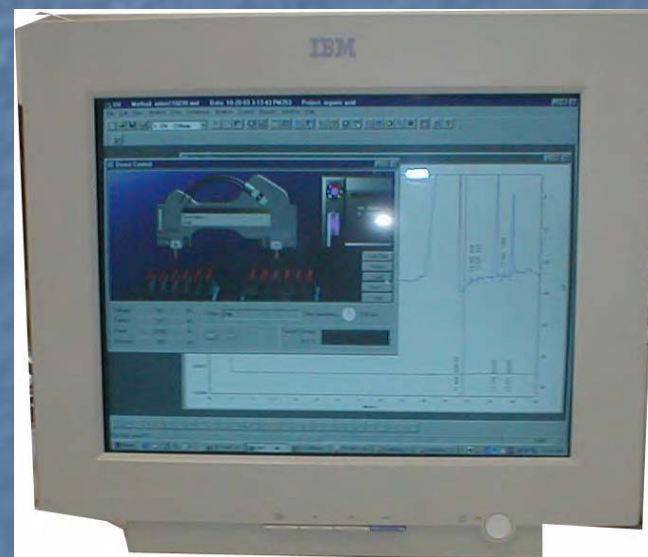



UV - 200nm: (0.00/15.00 Minutes) -- Current: (0.00/15.00 Minutes)

Time: 14.0014 Minutes - Amplitude: -- AU

Capillary

Detector

Samples and Buffers

25.0 °C
BI:B3 BO:B6

Description :
Lot# :

nm Absorb.
200 0.00000


Auto Zero
Home
Load
Inject...
Stop

Voltage: 0.0 kV
Current: 0.0 µA
Power: 0.000 W
Pressure: 20.0 psi

Status: Pause Time Remaining: 0.00 min

Sample Storage: 25.0 °C

Power Supply Pressure

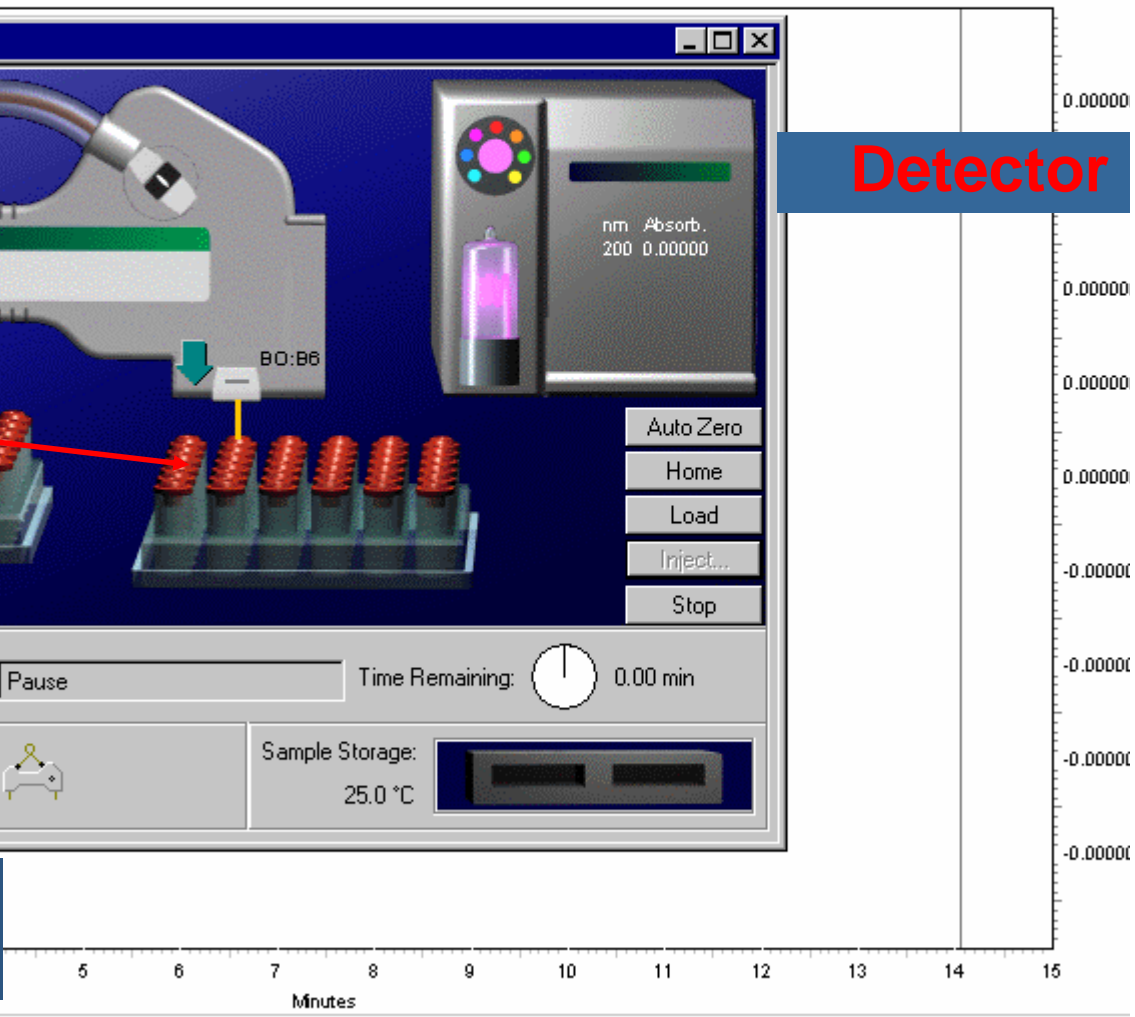




FRONT

BI




FRONT

BO

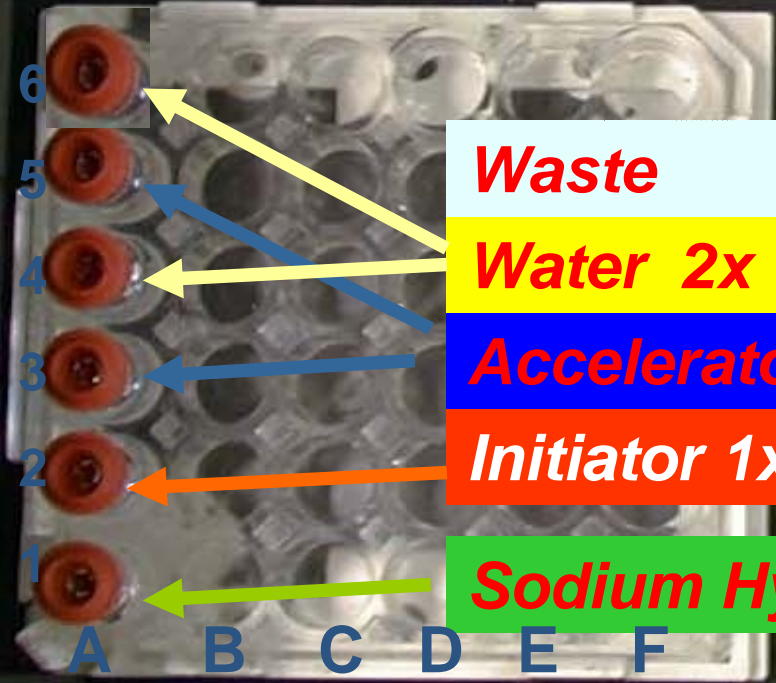

- Waste**
- Water 2x**
- Accelerator 3x 1.4mL**
- Initiator 1x 1.4mL**
- Sodium Hydroxide**

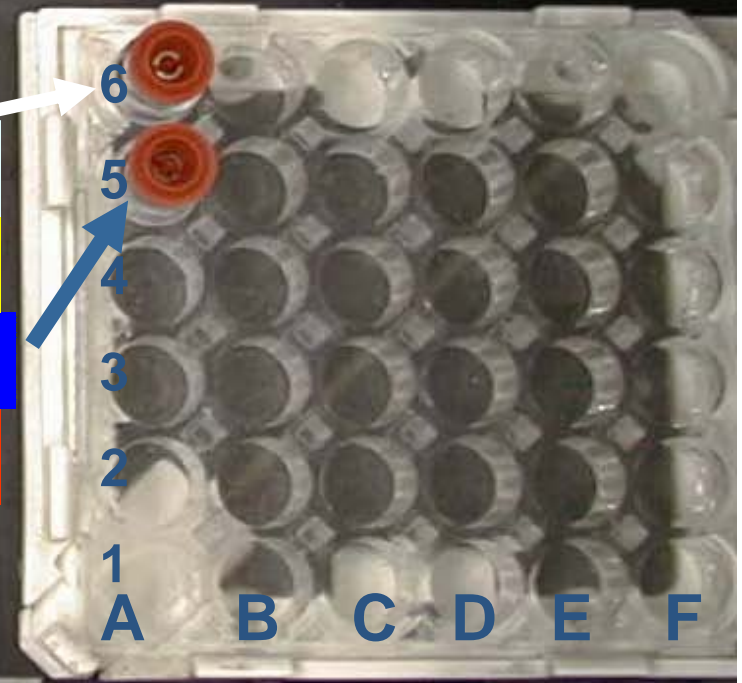


6
5
4
3
2
1
A B C D E F

CEofix™ Anions and Organic Acids


- Negative charge
 - CE: Reverse polarity
- Most are UV transparent
 - Buffer: indirect detection

CEofixTM Anions Organic Acids

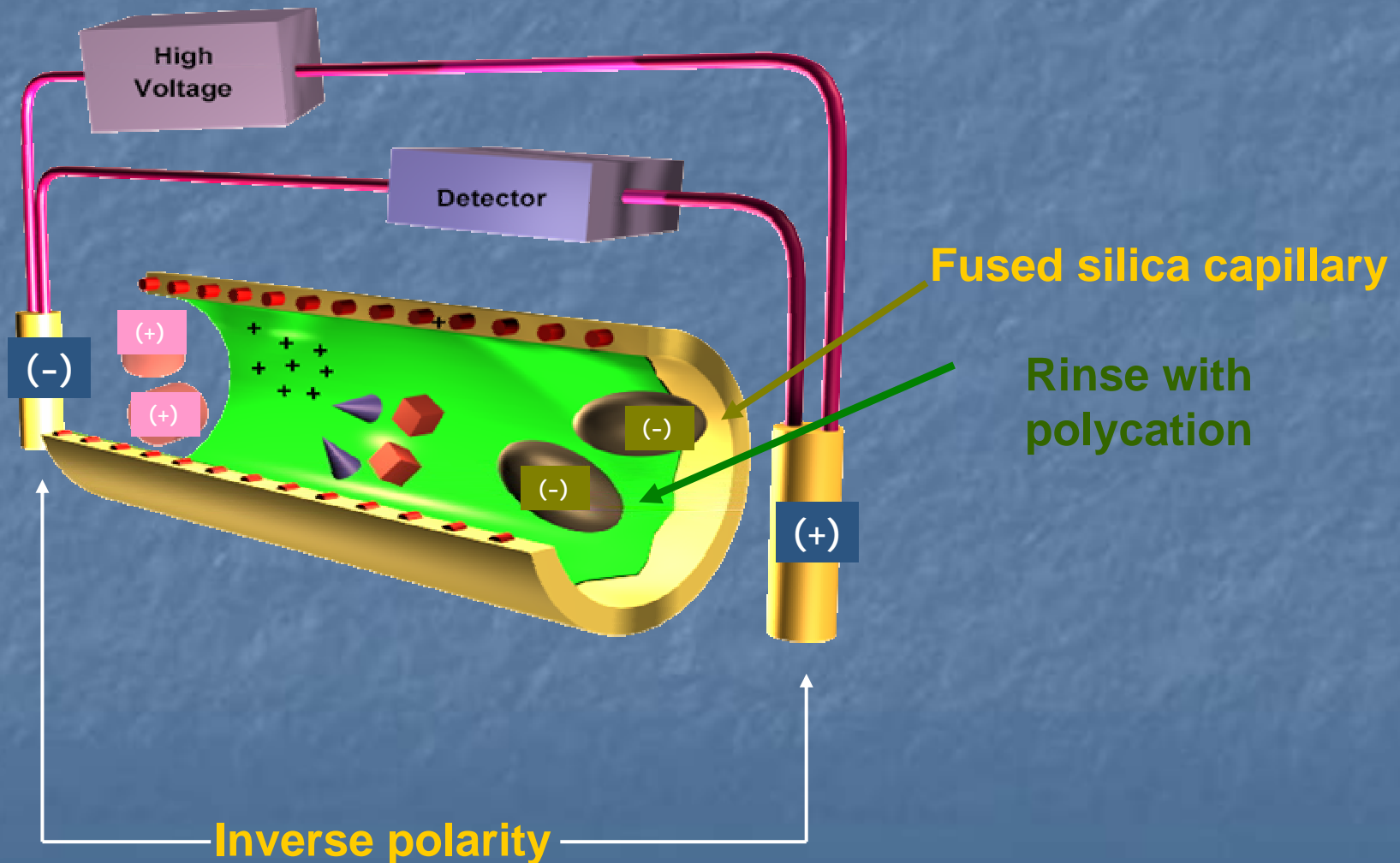

Detection: indirect-UV


P/ACE MDQ- DAD: 233 nm
UV filter at 230 nm
UV filter at 254 nm

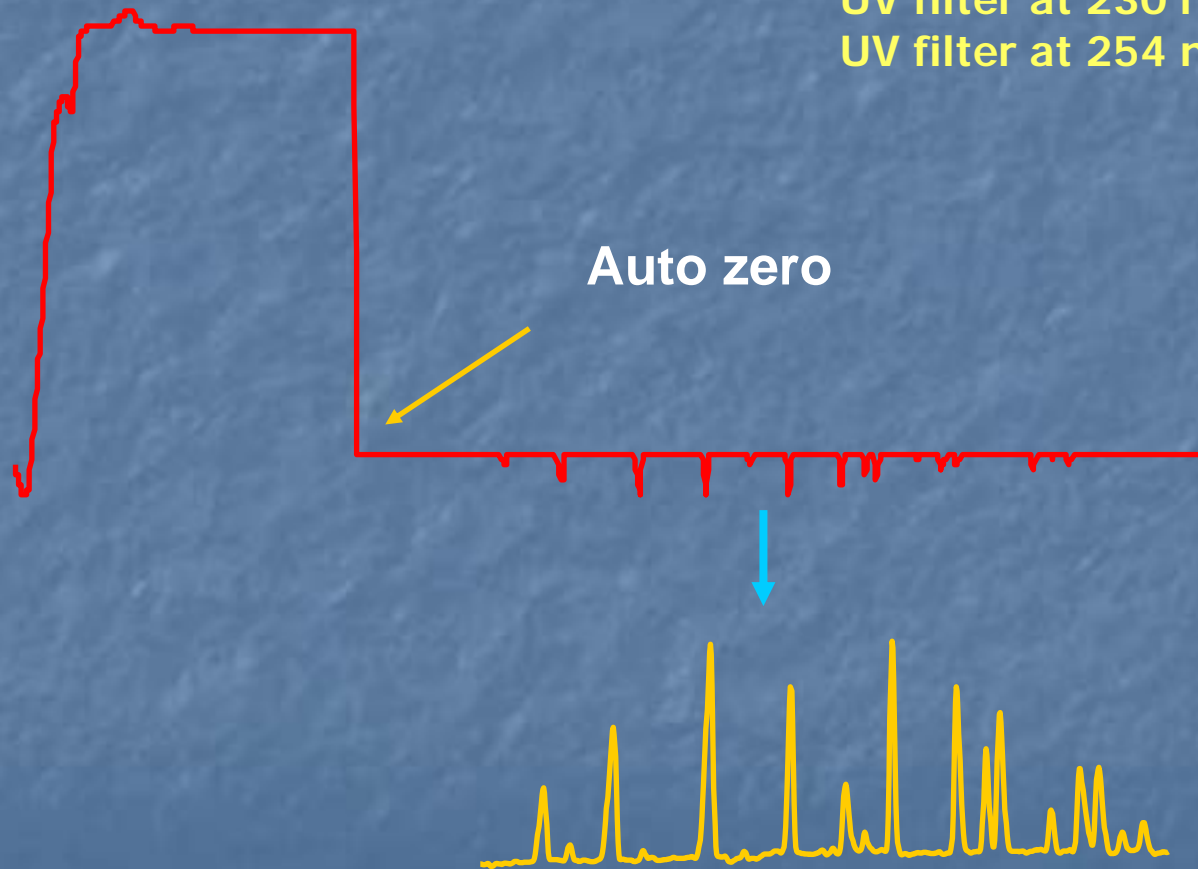

Composition


- **CEofix™ Anions5:**
pyridine-dicarboxylic pH 5.6
Reverse polarity, indirect detection (230 / 254 nm)
- **CEofix™ Anions8:**
pyridine-dicarboxylic pH 8.2
Reverse polarity, indirect detection (230 / 254 nm)
- **CEofix™ Anions2:**
phosphate pH 2.6
Reverse polarity and direct absorbance (200nm):
nitrate, nitrite, oxalate

Organic acids (CEofix Anions5) *pH 5.6*

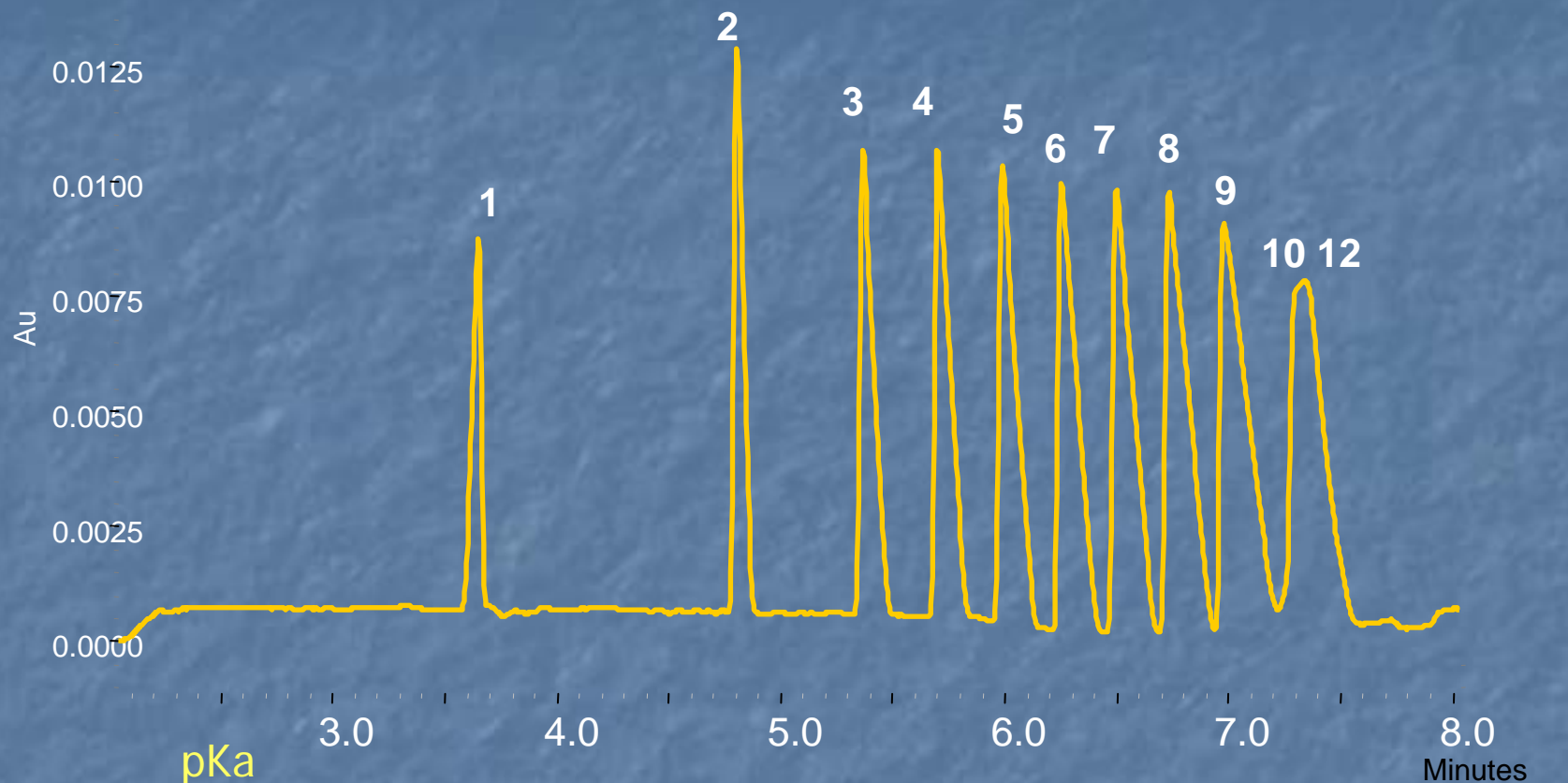

		pKa							
1	Formic	3.74	4	Butyric	4.87	7	Heptanoic	10	Capric
2	Acetic	4.76	5	Valeric	4.81	8	Caprylic	11	-
3	Propionic	4.87	6	Caproic	...	9	Pelargonic	12	Lauric

Organic acids

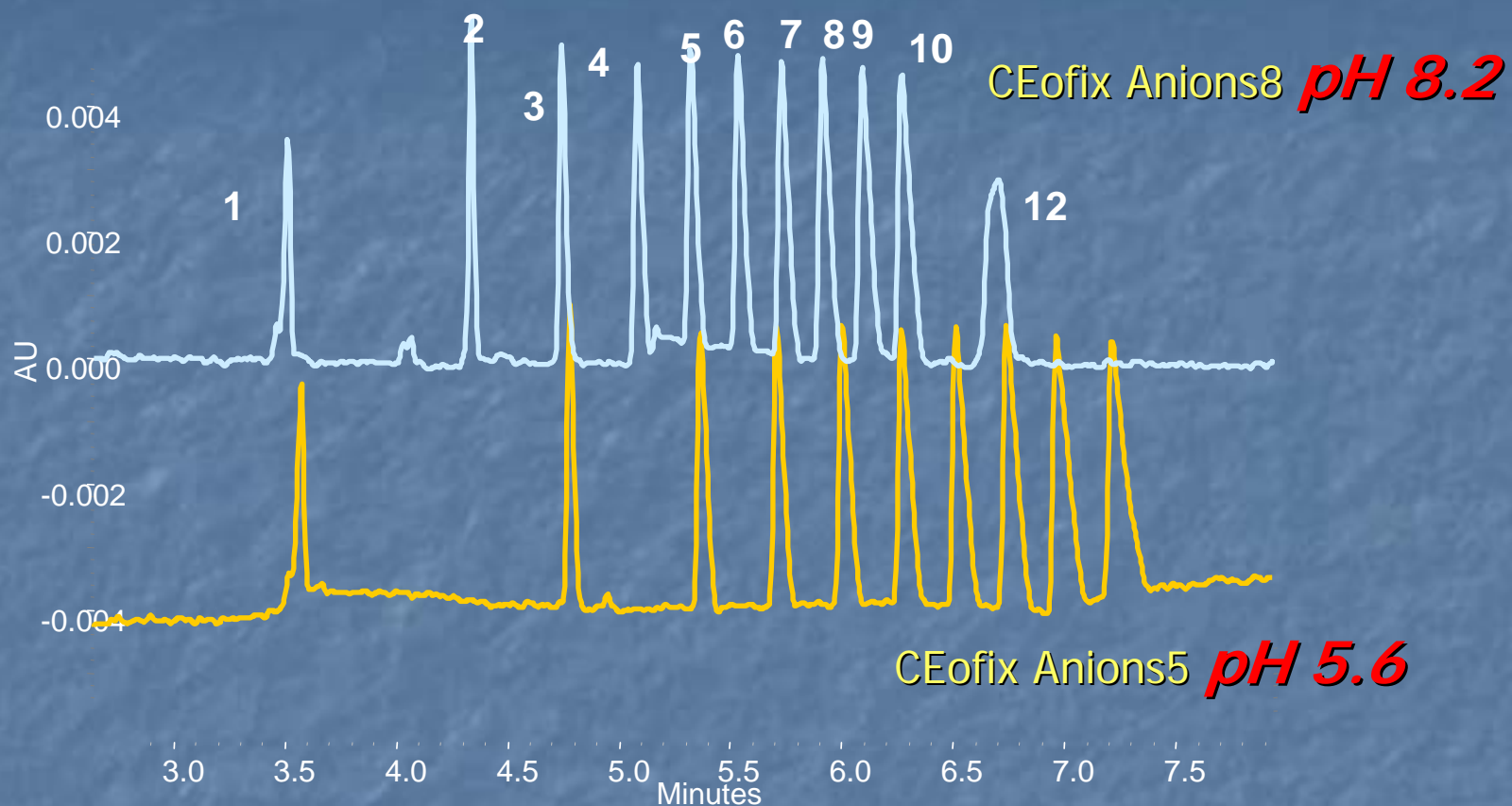

		pKa							
1	Formic	3.74	4	Butyric	4.87	7	Heptanoic	10	Capric
2	Acetic	4.76	5	Valeric	4.81	8	Caprylic	11	-
3	Propionic	4.87	6	Caproic	...	9	Pelargonic	12	Lauric

Test mix

- Anions:

- Chloride
- Nitrite
- Nitrate

- Organic Acid


- Propionic pKa 4.87
- Oxalic pKa: 1.25 – 3.67
- Malic pKa: 3.46 – 5.10
- Isocitric pKa: 3.29 – 4.71 – 6.40
- Citric: pKa: 3.13 – 4.76 – 6.40

Anions and Organic Acids: testmix


CEofix™ Anions 5
pyridine-dicarboxylic acid
(pH 5.6)



CEofix™ Anions 8
pyridine-dicarboxylic acid
(pH 8.2)

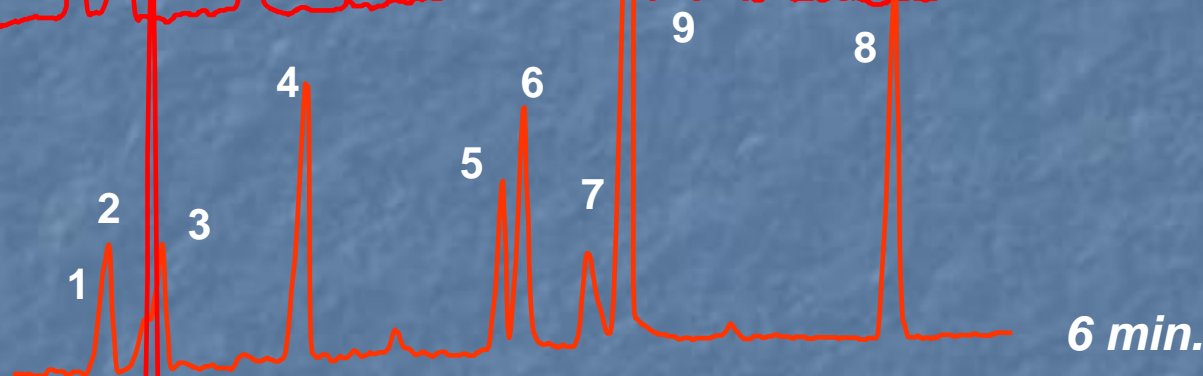

CEofix™ Anions 2
phosphoric acid
(pH 2.6)
direct UV


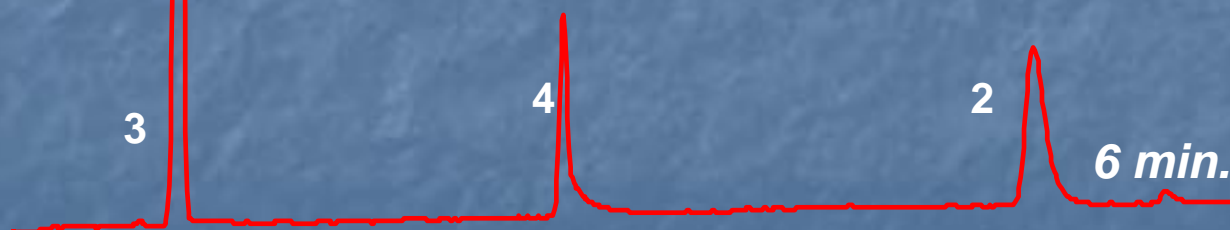


1. Chloride, 2. Nitrite, 3. Nitrate, 4. Oxalic, 5. Malic, 6. Iso-citric, 7. Citric, 8. Propionic, (9 Carbonate)

CEofix™ Cations


- Positive charge
 - CE: Normal polarity
- Some are UV transparent
 - Buffer: indirect detection
- CEofix™ Cation HR
 - 4-Aminopyridine pH 4.2*

CEofix™ Cations

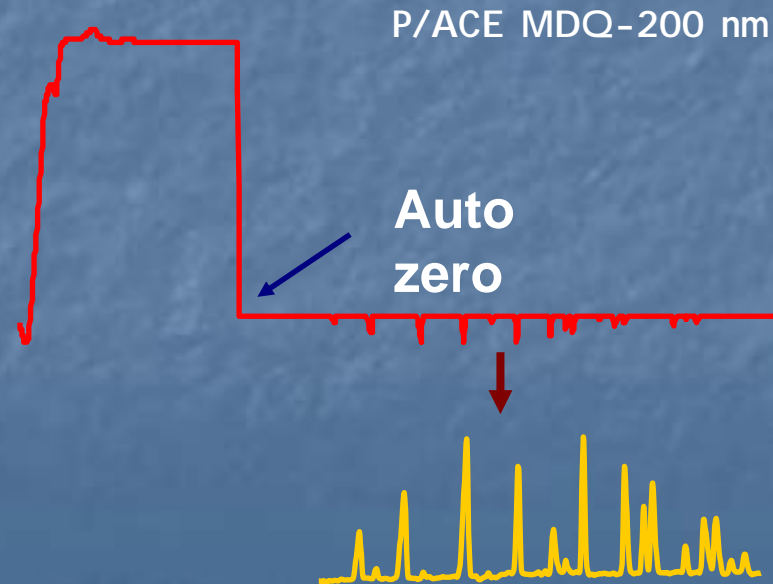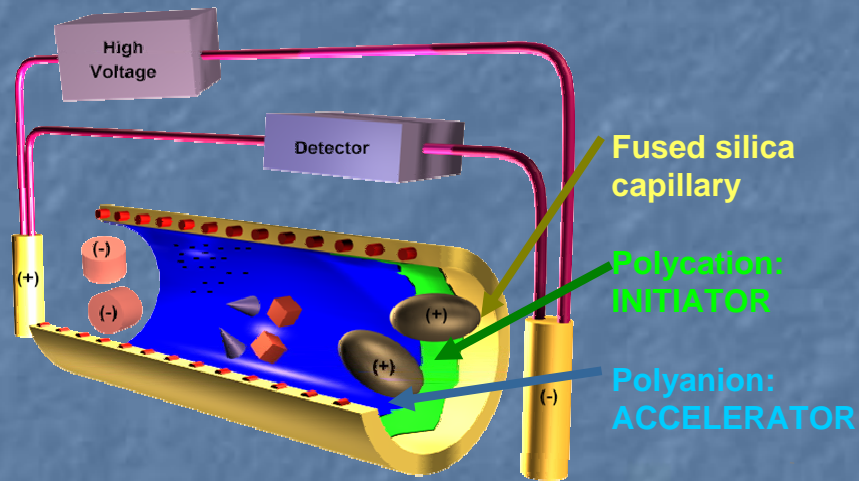

CEofix™ Cations


- Cations: 10 mg/L
(1) NH_4^+ , (2) K^+ , (3) Na^+ ,
(4) Li^+ , (5) Mg^{++} , (6) Ca^{++}
- Reproducibility on MT:
% RSD: < 0.5%
- Linearity: 1 to 25 mg/L
 $R^2 > 0.996$

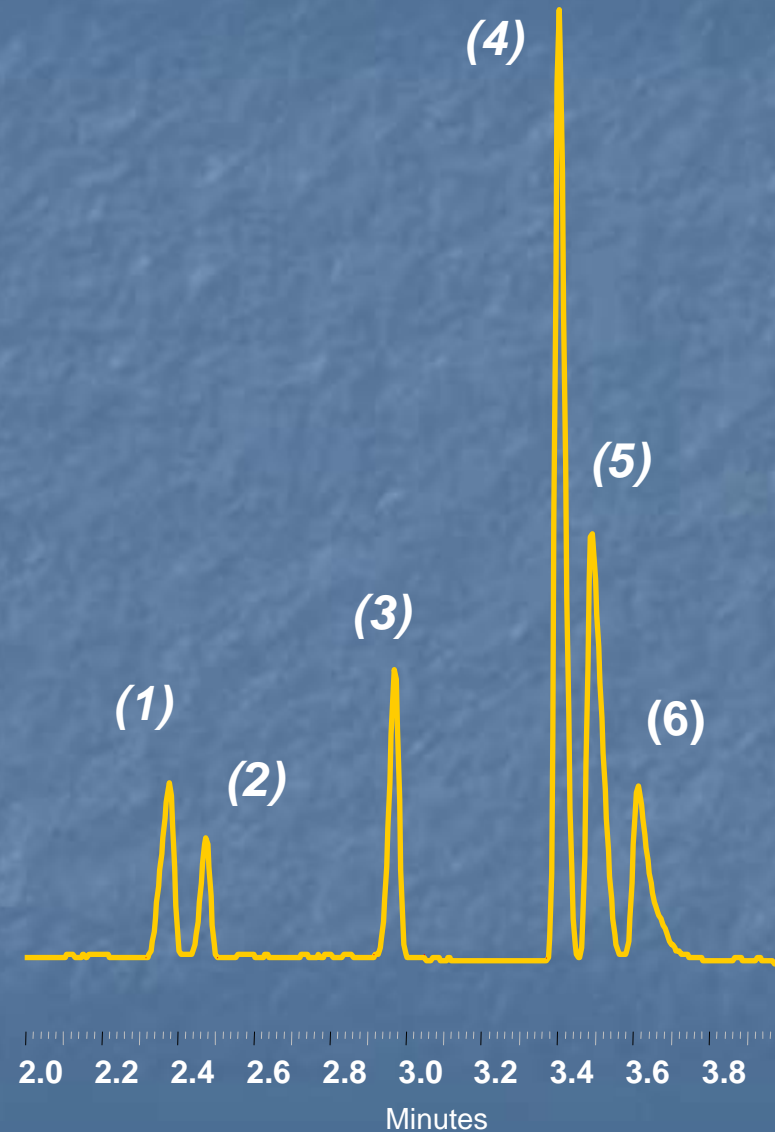

CEofix™ Cations

- Amines:

- (1) mono-ethanolamine,
- (2) di-ethanolamine, (3) N-benzyl-dimethylamine,
- (4) tri-ethanolamine, (5) 3-amino 1 phenylbutane,
- (6) dibenzylamine

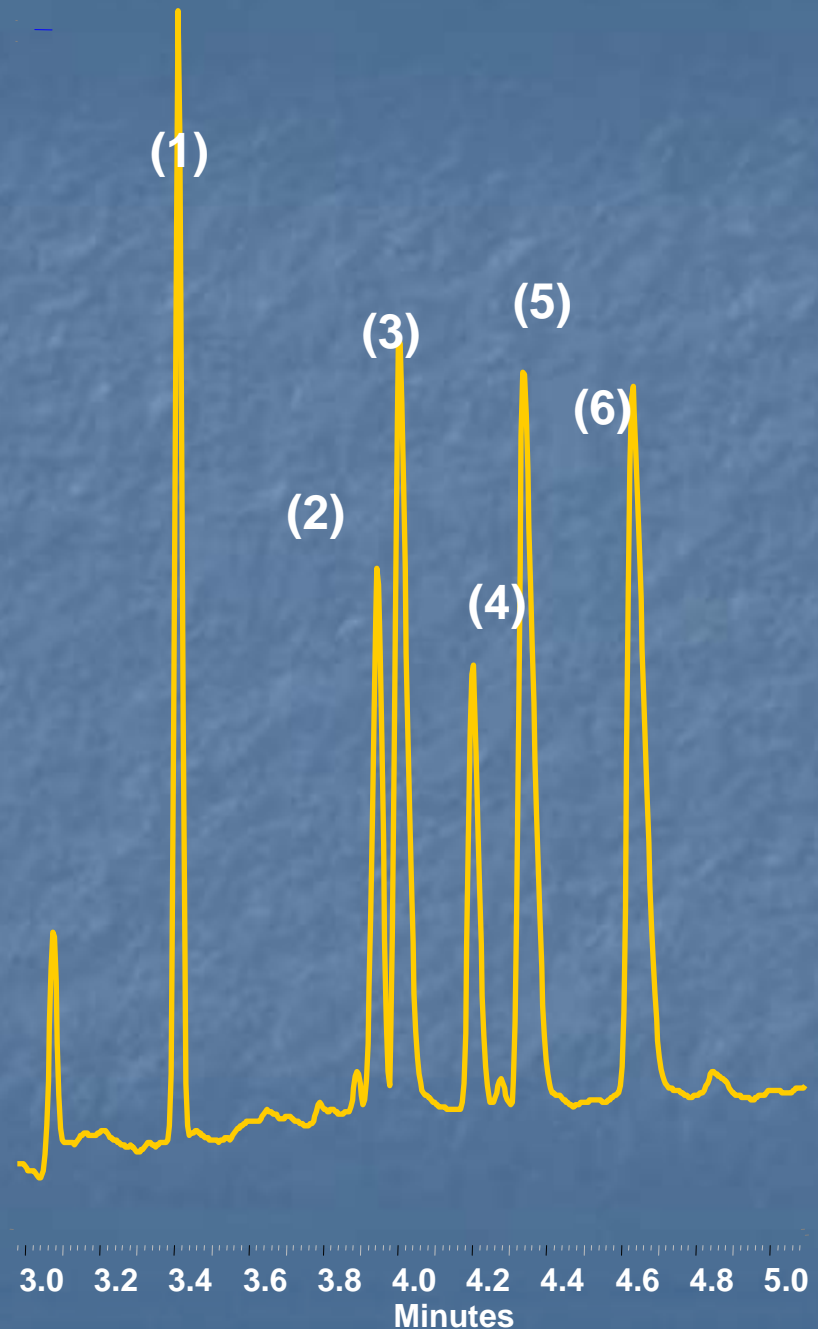


Applications

- General conditions
 - Beckman Coulter P/ACE MDQ
 - Capillary 75 μ x 60 cm (50cm to the detector)
 - Injection: 0.5psi during 5 sec.
 - Separation at 30 kV
 - Analysis time < 7 min
 - Detection UV filter or DAD

Application: urine

CEofix™ Anions5

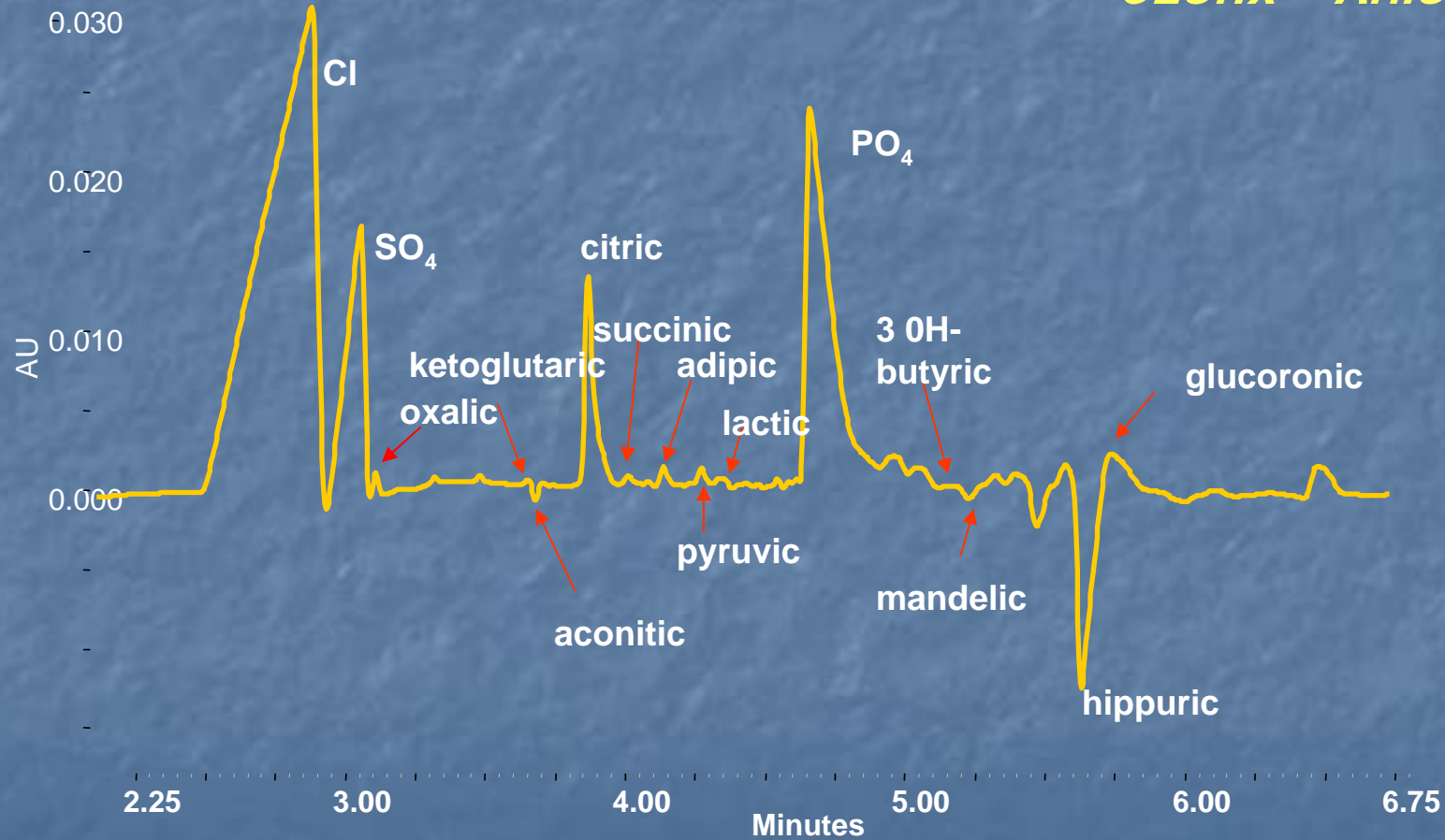

24 hrs - diluted ½ with H₂O

Application: Pharmaceutical Counter ions

- Drug: weak acid or base function
- Are produced as a salt
- Counter-ion quantification needed for P (purity)

$$P = 100\% - (m/M\% \text{ impurities} + m/M\% \text{ water} + m/M\% \text{ remaining solvents} + m/M\% \text{ inorganic material} + m/M\% \text{ Counter-ion})$$

For example: Sufentanyl citraat
total MW 578.69 \rightarrow 386.56 + 192.13
factor acid/base = $578.69/386.56 = 1.497$

Table 1

ORGANIC ANIONS MT in min.	Anions 5	Anions 2
Acetate	4.71	
Benzoate		10.36
Citrate	4.21	
Camphorsulfonate	5.89	
Di-benzoyl-tartrate		5.07
Di-p-Toluoyl-tartrate		5.38
Fumarate		6.71
2-Hydroxybenzoate		6.43
Hydroxybutanedioate	3.93	
2-Hydroxypropanoate	4.85	
Maleate		3.99
Malonate	3.89	
Mandelate		8.40
Methanesulfonate	3.99	
4-methylbenzenesulfonate		4.27
Oxalate	3.28	3.80
Propanoate	5.23	
Succinate	4.28	
Tartrate	3.73	

Table 2

INORGANIC ANIONS MT in min.	Anions 5	Anions 2
Nitrate HNO ₃	2.93	2.55
Phosphate H ₃ PO ₄	5.10	
Sulfate H ₂ SO ₄	3.11	
Chloride HCl	2.80	
Bromide HBr	2.77	2.40

Table 1 and 2 : Recorded migration time in minutes for the different anions, analysed with CEofix™ Anions5 and/or CEofix™ Anions2.

Table 3**Cations: Cations HR**


Ammonium (+1)

Calcium (+2)

Potassium (+1)

Magnesium (+2)

Sodium (+1)

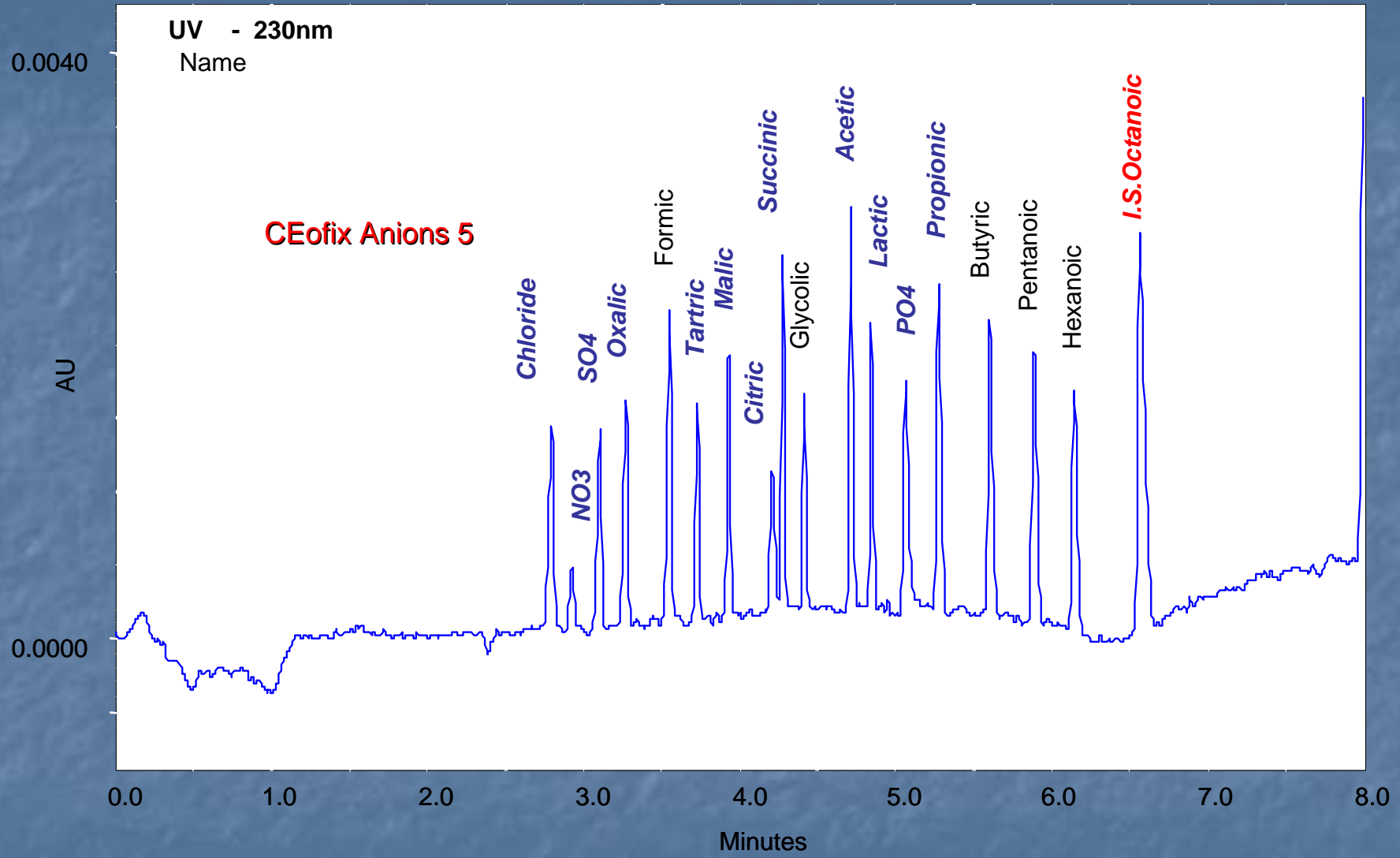

Succinic acid (-1) MT 4.13 / Succinate (-2) MT 4.11

Repeatability on migration time (octanoic acid as I.S.)

(N=10)	MT HCl	MT succinic acid	MT IS	Rel MT HCl	Rel MT succinic acid
	2.722	4.080	6.230	0.437	0.655
%SD	0.12%	0.07%	0.07%	0.09%	0.06%

Repeatability on area (octanoic acid as I.S.)


(N=10)	Area HCl	Area succinic acid	Area IS	Rel Area HCl	Rel Area succinic acid
	5523	3920	2218	2.490	1.768
%SD	2.91%	3.00%	3.22%	1.16%	1.06%

Silvie Verschakelen – Ivan Somers – Ivo Beyns – Rudy Sneyers – Achille Pluym

Johnson & Johnson Pharmaceutical Research & Development Drug Evaluation - Analytical Development

Application: animal feed

CEofix™ Anions5

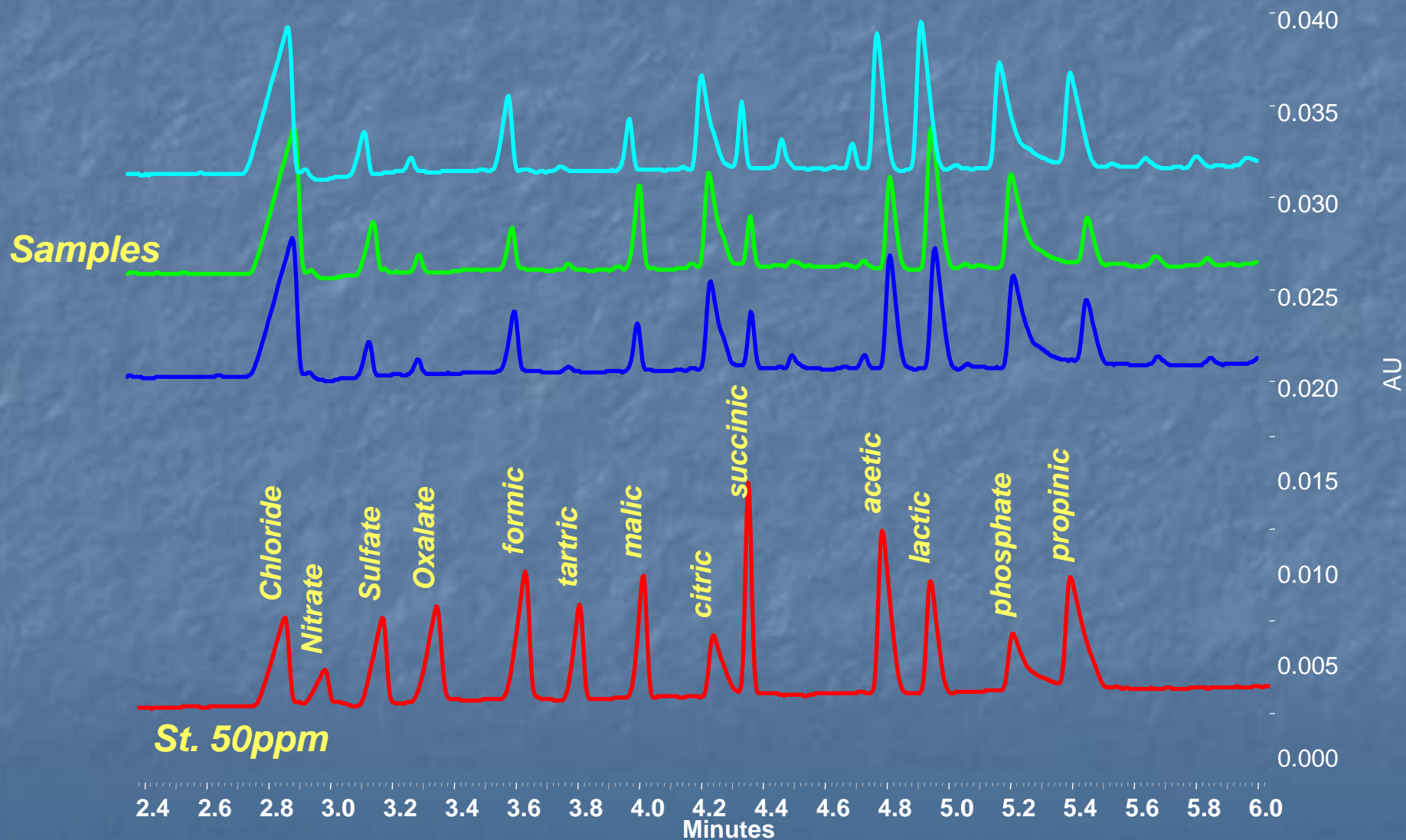


Application: Bromate in drinking water


CEofix Anions2 @ 200nm

Electrokinetic injection:
20 sec. @ 3 kV


R2 = 0.998

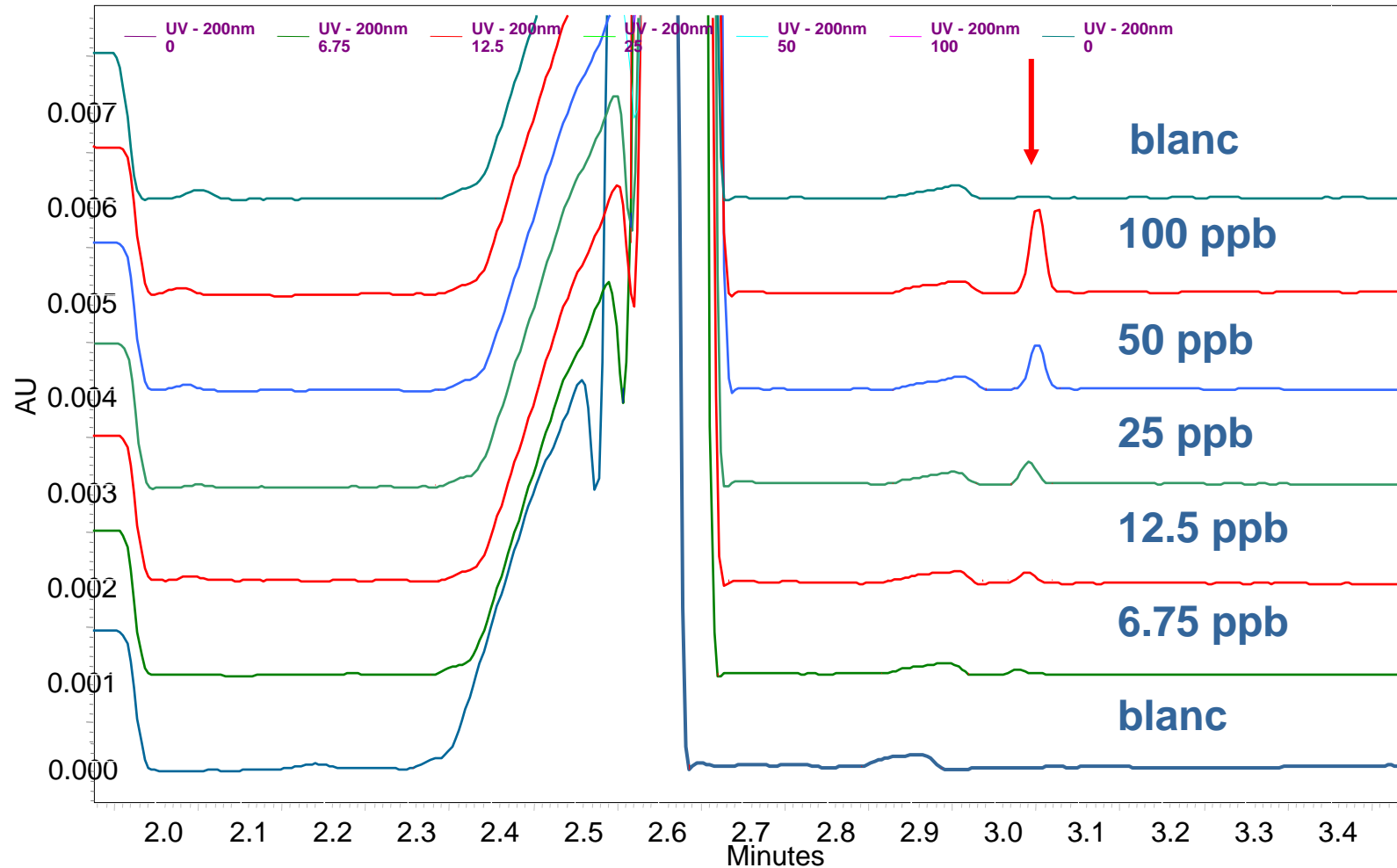

Application: mortar

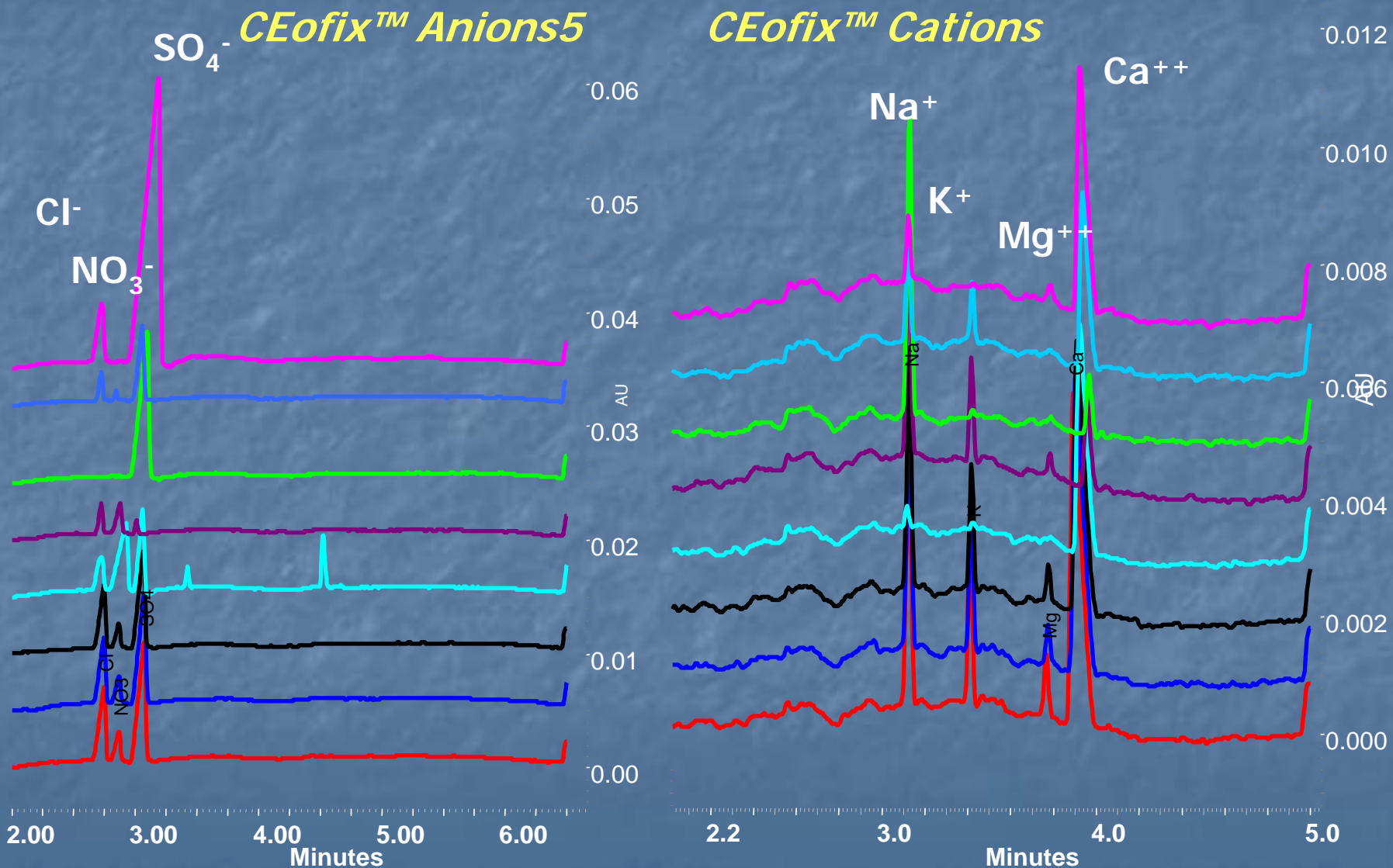

Applications: Amines in MeSO₃H (3.8 g/L)

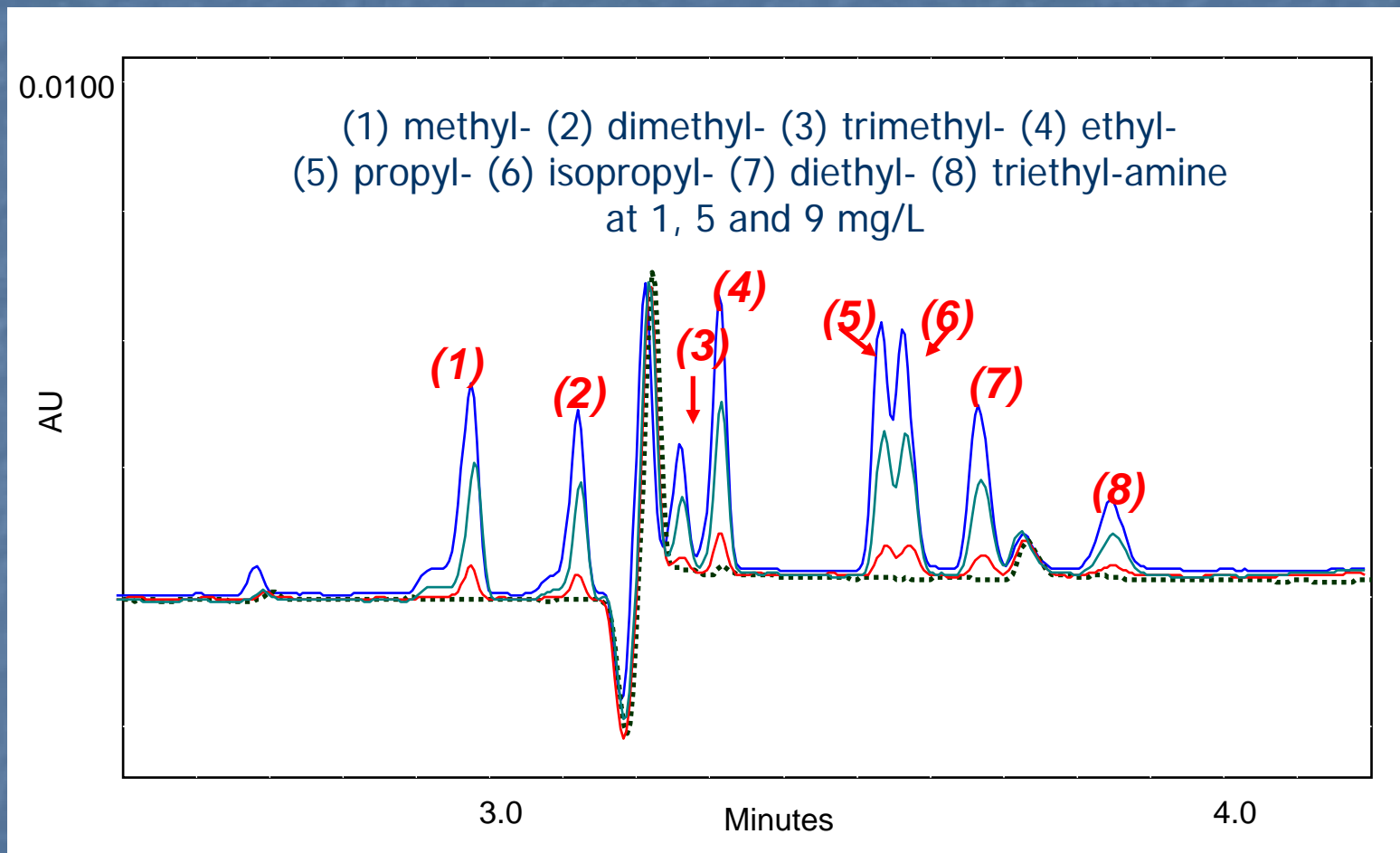

Buffer: diluted CEofix™ Cations HR – separation at 15°C

Points to consider when using CE

- No column but a capillary
- You work with ionised molecules
 - pH is very important
- Dilute your sample to be in the dynamic range
 - verify pH of dilution
- Be aware of matrix effects

Conclusion

- Fast method with CE
- Several applications on the same instrument
- CEofix™ buffers easy to use + support

ceofix@analisis.be

<http://www.analisis.be>

<http://ceofix.analisis.be/>